Ref. 100210



ES..... DESTILADOR / DESCALCIFICADOR.....	4
EN DISTILLER / DECALCIFIER.....	7
FR..... MACHINE À DISTILLER ET À DÉCALCIFIER.....	10
PO DESTILADOR / DESCALCIFICADOR.....	13
IT..... DISTILLATORE DECALCIFICATORE.....	16

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TALLERES MESTRAITUA S.L.
Txori-Erri Etorbidea, 60
Tfno. + 34 944530388 - Fax + 34 944711725
E-mail: mestra@mestra.es - www.mestra.es
48150 SONDIKA - BILBAO - ESPAÑA

Rev. 14/07/2022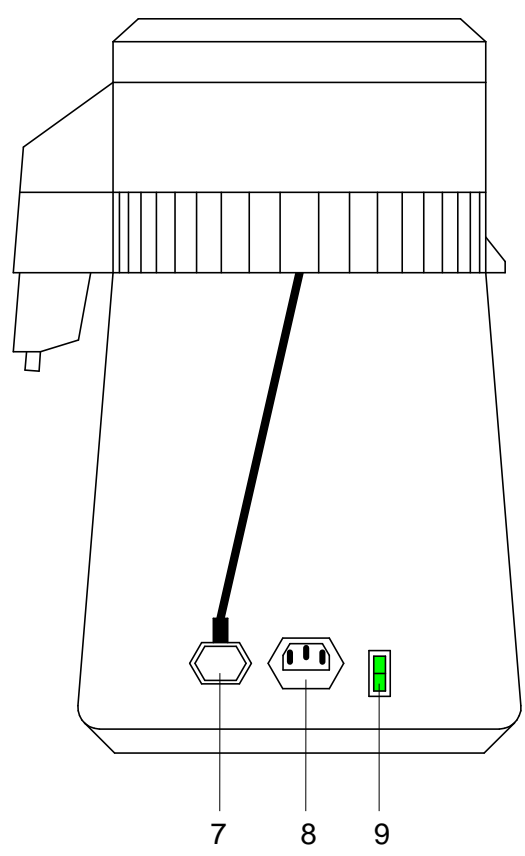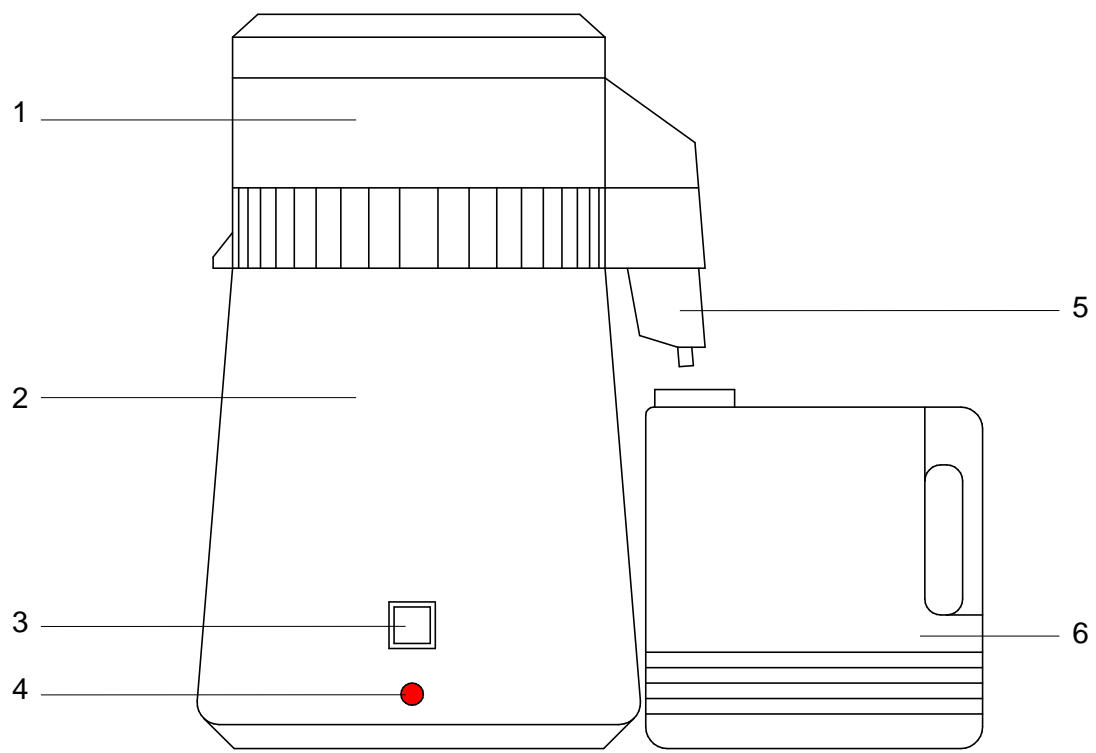

- 1 Tapa (ventilador) / **Lid (fan)** / *Couvercle (ventilateur)* / Tampa (Ventilador) / **Coperchio (ventilatore)**
- 2 Calderín / **Tank** / *Réservoir* / Reservatório / **Riscaldatore**
- 3 Reset
- 4 Piloto / **Pilot** / *Pilote* / Piloto / **Spia**
- 5 Salida agua / **Water outlet** / *Sortie d'eau* / Saída água / **Scarico acqua**
- 6 Garrafa / **Bottle** / *Goulot* / Boquilha / **Bocchello**
- 7 Cable ventilador / **Fan power cord** / *Câble ventilateur* / Fio ventilador / **Cavo ventilatore**
- 8 Conexión alimentación / **Power connection** / *Connexion alimentation* / Ligação alimentação / **Collegamento alimentazione**
- 9 Interruptor / **Switch** / *Interrupteur* / Interruptor / **Interruttore**

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The Ref. 100210 distiller/decalcifier is a simple and practical appliance: you will obtain 4 litres of distilled water by just filling the tank with tap water and waiting for about 6 hours. Its operation is fully automatic and silent. Cleaning the tank is easy and quick. The water produced is very pure (between 2 and 10 mg/litre of solid residue) and can be used in a large number of applications in the clinic and the dental lab (autoclaves, compound making, steam generating machines, etc.).

UNSUITABLE WATER FOR SURGICAL USE!

IMPORTANT- The appliance produces water of the "DISTILLED" type. It cannot be used in injections, vials, and wound cleaning or in any other surgical operation that requires "TRI-DISTILLED" type water.

SAFETY

- Read these instructions carefully before operating the unit.
- Keep these instructions for further consultation.

- Install the distiller on a dry, stable, heat-resistant and well-ventilated surface.
- The distiller requires a well-ventilated room. Bear in mind that while operating it emits heat.
- Use the unit exclusively for the aim it was designed.
- Do not allow children or non-qualified personnel handling the unit.
- Keep the citric acid container out of children's reach.

- Disconnect the unit from the power source for maintenance operations.
- Unplug the distiller for cleaning or refilling it.
- Do not dip any element of the distiller.
- If you need to interrupt the distillation cycle, stop the distiller by unplugging it.

- Do not open the distiller while it is working.
- Wait at least 20 minutes after a distillation cycle to open the distiller for cleaning it.
- Do not fill the tank with boiling water.

- **WARNING:** to prevent electric shock hazards, connect the unit to a proper power source (grounded 230 V, 50/60 Hz).

- Do not use cord extensions or connect any other appliance into the same outlet.
- Do not use the unit in damp rooms or near inflammable materials (alcohol, solvents...)

OPERATION

1. Put an activated carbon packet (L) in the nozzle. Rinse and clean it under the tap.

Note: fold the filter by half to place it in the cap.

2. Put the nozzle in its place.
3. Open the lid.
4. Pour water in the tank making sure not to exceed the mark. It is about 5 cm from the top.

Note: we recommend using hot water to shorten the time of the cycle.

5. Make sure all the contacts, wires and connections are not wet.
6. The lid of the bottle must be in line with the outlet hole of the tank. Plug the fan power cord and the main power cord in. Connect the main switch ON. You will notice that the fan starts working.

DO NOT ALLOW THE UNIT RUNS OUT OF WATER.

7. Turn off the distiller before the unit runs out of water in the kettle. This will prolong the life of the appliance.
8. However, the distiller will automatically shut off after the cycle is complete. In that case, press the “RESET” button to restart the distillation.

CLEANING THE TANK

- Fill the tank with 4 litres of hot water.
- Add 1-3 spoonfuls of cleaner (too much acid may damage the material of the tank).
- Leave the lid open, connect the power cord in and press the switch to heat the tank until the water boils.
- Empty the tank after 15-20 minutes. Then, wash the inside with clean water.
- If there is too much residue to be removed in one go, repeat the process several times until the inside is completely clean.
- A small amount of residue can cause no damage and you can use the unit safely.

DO NOT ADD ACID TO WATER. The acid will not evaporate, will increase the acidity cycle after cycle and the it will finally perforate the tank.

MAINTENANCE

- **Do not allow the unit runs out of water.** Turn off the distiller before the unit runs out of water in the kettle. This will prolong the life of the appliance.
- **DO NOT ADD ACID TO WATER.**
- Do not use abrasive cleaners or brushes. Use only a specific cleaner for distillers.
- Clean the tank approximately once a week or more often if the water is hard.
- Allow enough time for the cleaner to work and dissolve the particles set into the tank.
- The filter pack should be replaced at least once a month.
- If the coil is dirty, the distillation process will be less efficient. Clean the coil regularly through the upper grille with a brush and suction pump.
- After each operation, rinse the tank to prevent the residues hardening.
- Keep the outside of the unit dry.

TECHNICAL SPECIFICATIONS

Voltage:	230 V, 50/60 Hz
Power:	580 W
Resistance:	565 W
Fan:	15 W
Distillating capacity:	approx. 4 L in 6 h
Height:	380 mm
Width:	250 mm
Depth:	260 mm
Tank capacity:	4 L
Weight:	3,5 kg

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